Based on the principle of positive displacement, StepMate Stepper is designed for trouble-free and reliable dispensing of liquids, especially for high viscosity, high density and high vapor pressure. The 48-step dispensing allows for enhanced accuracy and usability.

StepMate

Stepper

Features

- Light and ergonomic solution provides easy usability
- Single-handed operation
- Volume range from 10µL to 5mL
- Up to 48 dispensing steps
- Maintenance free
- Equipped with durable tip
- insertion lever
- Fits seven sizes of disposable polypropylene syringes from 0.5mL to 50mL

How to select dispensing volume

Find the required dispensing volume in the table Use the adjusting wheel to set the required volume Select and insert a suitable tip Follow the maximum dispensing steps

Tip volume

Di	Dispensing Volume									
Tips Sel.	1	2	3	4	5					
0.50ml	10µl	20µI	30µl	40µI	50µl					
1.25ml	25µI	50µI	75µI	100µI	125µI					
2.50ml	50µl	100µl	150µl	200µl	250µI					
5.00mi	100µl	200µI	300µl	400µl	500µl					
12.50ml	250µl	500µl	750µl	1000µl	1250µl					
25.00ml	500µI	1000µl	1500µl	2000µl	2500µl					
50.00ml	1000µl	2000µl	3000µl	4000µi	5000µl					
Steps	48	23	15	11	8					

Steps



Optimized for smooth dispensing of even the smallest volumes

Five easy-to-select settings control

dispensing volume



20

tepMate

Stepper

DLAB stepper tips are particularly suitable for dispensing of fluids with high viscosity, high vapor pressure etc.

Tips are made from high-quality materials (Cylinder: PP, Piston: PE-HD) without soft lubricants or other additives.

1 2 3 4 5

Specifications

Dial setting

Note: The first and last dispensing liquid should be discarded for precision.

<u></u>			! *	! *	! +	: *	4	2 Martin	
Number o	f steps	48	23	15	11	8			
Cat. No.	Syringe volume (mL)	Sample volume (µL)					Α%	Р%	Description
17900032	0.50	10	20	30	40	50	± 0.8	≤ 0.7	0.5mL, piston PE-HD, cylinder PP, 1 packing, 5 pcs/packing
17900033	1.25	25	50	75	100	125	± 0.8	≤ 0.5	1.25mL, piston PE-HD, cylinder PP, 1 packing, 5 pcs/packing
17900034	2.50	50	100	150	200	250	± 0.8	≤ 0.4	2.5mL, piston PE-HD, cylinder PP, 1 packing, 5 pcs/packing
17900035	5.00	100	200	300	400	500	± 0.4	≤ 0.3	5mL, piston PE-HD, cylinder PP, 1 packing, 5 pcs/packing
17900036	12.50	250	500	750	1000	1250	± 0.3	≤ 0.3	12.5ml, piston PE-HD, cylinder PP, 1 packing, 5 pcs/packing
17900037	25.00	500	1000	1500	2000	2500	± 0.2	≤ 0.3	25mL, piston PE-HD, cylinder PP, 1 packing, 2pcs/packing with 1 pc adapter
17900038	50.00	1000	2000	3000	4000	5000	± 0.2	≤ 0.2	50mL, piston PE-HD, cylinder PP, 1 packing, 2pcs/packing with 1 pc adapter

Stepper

2020/11/30 上午10:36